**YS Series
Instruction Manual**

**Serie YS
Manual de instrucciones**

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1. INTRODUCTION

This manual contains installation, operation and maintenance instructions for the YS Series Scales. Please read the manual completely before using the scale.

1.1 Safety Precautions

Please follow these safety precautions:

- Verify that the AC Adapter input voltage matches the local AC power supply.
- Use the scale only in dry locations.
- Do not operate the scale in hostile environments.
- Do not drop loads on the platform.
- Do not place the scale upside down on the pan.
- Service should be performed only by authorized personnel.

1.2. Controls

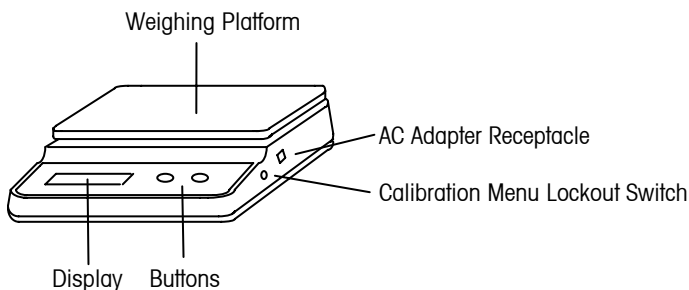
Figure 1-1. Controls.

TABLE 1-1. CONTROL FUNCTIONS.

Button	Functions
TARE <i>On</i> Yes	Short Press: (when Off) Turns scale on
	(when on) Sets display to zero
	Tares weight of item on pan
MODE <i>Off</i> No	Long Press: Enters menu
	Short Press (Menu) Selects the displayed item
MODE <i>Off</i> No	Short Press Steps through active units and modes
	Long Press: Turns the scale Off
	Short Press (Menu) Advances to the next selection

2. INSTALLATION

2.1 Package Contents

- Scale
- Pan
- AC Power Adapter
- Instruction Manual
- Warranty Card

2.2. Location

Use the scale on a firm, steady surface. Avoid locations with excessive air current, vibrations, heat sources, or rapid temperature changes.

2.4 Power

Power the scale using the AC power adapter (included) or 4 AA batteries.

AC Power Connection

Plug the jack from the adapter into the receptacle located on the right side of the balance. Connect the AC adaptor to a properly grounded electrical outlet.

Battery Installation

Turn the balance over and remove the battery cover. Install 4 AA batteries in the battery compartment, orienting the minus (-) ends against the coil springs and the plus (+) ends against the buttons. Replace the battery cover.

2.5 Initial Calibration

When the scale is first installed, it should be calibrated to ensure accurate weighing results. Refer to section 4.1 Calibration Menu.

3. OPERATION

3.1 Weighing Mode

Use this mode to measure the weight of a sample in the selected unit of measure.

1. Repeatedly press **MODE** until the desired unit is displayed.
2. Press **TARE** to set the display to zero.
3. Add material to the pan.
4. The display shows the weight of the material on the pan.

To weigh material in a container:

1. Press **TARE** to set the display to zero.
2. Put an empty container on the pan; then press **TARE** to tare the container weight.
3. Add material to the container.
4. The display shows the net weight of the material.

3.2 Counting Mode

Use this mode to count parts of uniform weight. The balance determines the quantity based on the average weight of a single part. All parts must be uniform in weight for accurate measurements.

1. Repeatedly press **MODE** until PC **CON** is displayed.
2. Put an empty container on the pan; then press **TARE** to tare the container weight.
3. The display shows the sample size PC **Add 5** .
4. To change the sample size, repeatedly press **TARE/On** until the desired quantity is displayed: PC **Add 10**, PC **Add 20**, PC **Add 30**, PC **Add 40**, PC **Add 50**, PC **Add 5...**
5. Add the displayed number of parts to the container; then press **TARE/On**.
6. The display shows the current number of parts PC **5** .

Note: To exit or restart Counting at any time, press and hold **TARE/On** until PC **CON** is displayed. Return to step 2 to restart, or press **MODE/Off** for other weighing functions.

4. MENUS

The menu allows the user to calibrate the scale and to customize scale settings.

To enter the menu, press and hold **TARE/On/Yes** button. To enter the Calibration menu, release the button when **CAL** is displayed. To enter the Setup menu, continue holding the button and release it when **SETUP / RESET** is displayed.

Note: The Calibration Lockout Switch must be in the "UNLOCK" position to enter the Calibration menu.

4.1 Calibration Menu [CAL]

Upon entering the Calibration menu, the Calibration Weight setting is displayed.

Press **Yes** to accept the setting or press **No** to change the setting.

TABLE 2-1. CALIBRATION WEIGHTS.

Model	YS202	YS302	YS601	YS1101	YS2101	YS5001
Default Calibration Weight	200 g	300 g	600 g	1000 g	2000 g	5000 g
Alternate Calibration Weight	100 g	100 g	300 g	N/A	1000 g	N/A

When the desired Calibration Weight setting is displayed, press **Yes**. The display flashes 0, momentarily. Then the display flashes the Calibration Weight value. Place the Calibration Weight on the pan. When the display stops flashing, calibration is completed and the scale returns to the weighing mode.

Note: To prevent unauthorized calibration, position the Calibration Lockout Switch in the "LOCK" position.

4.2 Setup Menu [SETUP]

To navigate through the Setup menu, press **Yes** to view the setting for the menu item or press **No** to move to the next menu item. When viewing the setting, press **Yes** to accept the setting, or press **No** to change the setting.

RESET [RESET]

Reset the Setup menu to the factory defaults.

no = not reset

YES = reset

UNIT SELECTION [SEL]

Set the status of each weighing unit.

OFF = unit is disabled

ON = unit is enabled

The symbol for each unit is displayed next to the setting (g = grams, oz = ounces, ozt = troy ounces, dwt = pennyweights, lb = pounds, t = taels, ct = carat, PC = pieces).

When t is set to \square , select the type of taels (t = Hong Kong, t = Singapore or t = Taiwan).

STABILITY RANGE [StAb]

Set the stability range.

- .5 d = the unit symbol will light when the display is stable within 0.5 divisions
- 1 d = the unit symbol will light when the display is stable within 1 division
- 2 d = the unit symbol will light when the display is stable within 2 divisions
- 5 d = the unit symbol will light when the display is stable within 5 divisions

AUTO ZERO TRACKING RANGE [AZTR]

Set the automatic zero tracking range.

- .5 d = the display will maintain zero until a drift of 0.5 divisions per second has been exceeded
- 1 d = the display will maintain zero until a drift of 1 division per second has been exceeded
- 3 d = the display will maintain zero until a drift of 3 divisions per second has been exceeded
- OFF = disabled

PARTS COUNTING ERROR LEVEL [PCER]

Set the parts counting error level.

- OFF = PCER message is disabled
- 1 PCt = PCER is displayed if the count accuracy exceeds 1 percent
- 2.5 PCt = PCER is displayed if the count accuracy exceeds 2.5 percent
- 5 PCt = PCER is displayed if the count accuracy exceeds 5 percent

AUTO OFF [A. OFF]

Set the automatic shut off timer.

- OFF = disabled
- 60 = the scale turns off after 60 seconds of inactivity
- 120 = the scale turns off after 120 seconds of inactivity
- 180 = the scale turns off after 180 seconds of inactivity

END [End]

Press **Yes** to exit the Setup menu or **No** to return to the top of the Setup menu.

5. MAINTENANCE

5.1 Troubleshooting

The following table lists common problems and possible causes and remedies. If the problem persists, contact Ohaus or your authorized dealer.

TABLE 5-1. TROUBLESHOOTING.

Symptom	Possible Cause	Remedy
Cannot turn on	No power to scale	Verify connections and voltage.
Poor accuracy	Improper calibration • Unstable environment	Perform calibration • Move scale to suitable location
Cannot calibrate	Unstable environment • Incorrect calibration weight	Move the scale to suitable location • Use correct calibration weight
Cannot access unit	Unit not enabled	Enter menu and enable unit
Err 0	Pan was removed during weighing	Re-install pan
Err 1	Weight on pan exceeds capacity	Remove weight from the pan
Err 2	Pan was removed prior to power on	Install pan and re-zero.
Err 3	Pan has load during power on	Remove weight from pan and re-zero.
PCErr	Counting sample size is too small	Use larger sample
Low battery indicator	Batteries are weak	Replace batteries

5.2 Service Information

If the troubleshooting section does not resolve or describe your problem, contact your authorized Ohaus service agent. For service assistance or technical support in the United States call toll-free 1-800-526-0659 between 8:00 AM and 5:00 PM EST. An Ohaus product service specialist will be available to provide assistance. Outside the USA, please visit our web site, www.ohaus.com to locate the Ohaus Office nearest you.

6. TECHNICAL DATA

The technical data is valid under the following ambient conditions:

Ambient temperature: 10°C to 40°C

Relative humidity: Maximum 80% up to 31°C, decreasing linearly to 50% at 40°C

Operability is assured at ambient temperatures between 10°C and 40°C

Power: AC adapter 6 VDC 200 mA output

Protection: Protected against dust and water

Pollution degree: 2

Installation category: Class III

EMC: See Declaration of Conformity

6.1 Specifications

TABLE 6-1.SPECIFICATIONS.

Model	YS202	YS302	YS601	YS1101	YS2101	YS5001
Capacity (g)	200	300	600	1100	2100	5000
Readability (g)	0.01	0.05	0.1			0.5
Repeatability (g)	0.015	0.075	0.15			0.75
Linearity (g) (+/-)	0.02	0.1	0.2			1
Weighing Units	grams, ounces, troy ounces, pennyweights, pounds, taels (3), carats					
Application Modes	Weighing, Counting					
Tare range	To capacity by subtraction					
Stabilization Time	3 seconds					
Power Requirements	6 VDC 200 mA AC adaptor (supplied) or 4 AA (LR6) batteries (not supplied)					
Calibration	Digital with external weight					
Display Type	5-digit 7-segment LCD					
Display Size (mm / in)	12.7 / 0.5 digits					
Keypad	2-button overlay					
Overall Size (mm / in)	176 x 60 x 194 / 6.9 x 2.4 x 7.6					
Pan Size (mm / in)	120 / 4.7		166 x 128 / 6.5 x 5			
Net Weight (kg / lb)	0.8 / 1.8					
Shipping Weight(kg/lb)	1.5 / 3.3					

6.2 Compliance

Compliance to the following standards is indicated by the corresponding mark on the product.

Mark	Standard
	This product conforms to the EMC directive 2004/108/EC and the Low Voltage Directive 2006/95/EC. The complete declaration of Conformity is available from Ohaus Corporation.

	<p>Disposal</p> <p>In conformance with the European Directive 2002/96 EC on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.</p> <p>Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment.</p> <p>If you have any questions, please contact the responsible authority or the distributor from which you purchased this device.</p> <p>Should this device be passed on to other parties (for private or professional use), the content of this regulation must also be related.</p> <p>Thank you for your contribution to environmental protection.</p> <p>For disposal instructions in Europe, refer to www.ohaus.com/weee</p>
---	---

FCC Note

This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.

Industry Canada Note

This Class B digital apparatus complies with Canadian ICES-003.

ISO 9001 Registration

In 1994, Ohaus Corporation, USA, was awarded a certificate of registration to ISO 9001 by Bureau Veritas Quality International (BVQI), confirming that the Ohaus quality management system is compliant with the ISO 9001 standard's requirements. On May 15, 2003, Ohaus Corporation, USA, was re-registered to the ISO 9001:2000 standard.

LIMITED WARRANTY

Ohaus products are warranted against defects in materials and workmanship from the date of delivery through the duration of the warranty period. During the warranty period Ohaus will repair, or, at its option, replace any component(s) that proves to be defective at no charge, provided that the product is returned, freight prepaid, to Ohaus. This warranty does not apply if the product has been damaged by accident or misuse, exposed to radioactive or corrosive materials, has foreign material penetrating to the inside of the product, or as a result of service or modification by other than Ohaus. In lieu of a properly returned warranty registration card, the warranty period shall begin on the date of shipment to the authorized dealer. No other express or implied warranty is given by Ohaus Corporation. Ohaus Corporation shall not be liable for any consequential damages.

As warranty legislation differs from state to state and country to country, please contact Ohaus or your local Ohaus dealer for further details.

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P/N 80252149 © 2008 Ohaus Corporation, all rights reserved / todos los derechos reservados / tous droits réservés / Alle Rechte vorbehalten / tutti i diritti riservati.